

Phycocyanobilin (PCB) Purification from *Spirulina*

Obtain cyanobacteria, *Spirulina* (Seltzer Chemical(Spirulina Powder , 191775-C, 1-800-735-8137)

(<http://www.seltzerchemicals.com/Products/All/>)

Day 1

1 –Suspend 50 g *Spirulina* powder in 1.5 liter H₂O (30 ml/g) in a 2.8-liter flask covered with aluminum foil. Stir for 10 minutes then divide into six 250 ml centrifuge bottles and centrifuge 8000 RPM, GS3 rotor, 4° C for 1 hr.

2 – Decant supernatant into another 2.8 liter flask and add 15g TCA - final concentration of 1% (w/v). Stir for 1 hr in the dark (aluminum foil cover) at 4° C. Centrifuge 8000 RPM, GS3 rotor, 4° C for 10 min. Discard supernatant.

3 - Add 25 ml MeOH to each bottle; resuspend green pellets completely using a tissue homogenizer and/or vortex – keep pellets on ice during this process. After pellets are fully resuspended, add MeOH to each bottle to a final volume of 250 ml and repeat centrifugation (8000 RPM, 4° C / 10 min.) Discard supernatant.

4 – Repeat step (3) 4X – or until MeOH wash is colorless.

5 – Store final washed pellets at -20° C, wrapped in foil overnight.


******It is critical that all steps from this point on are performed under low intensity, green ‘safelight’ conditions in a darkroom (see below for safelight specifications - a photographic darkroom, with “orange” safelights turned off, is fine); and/or that samples are shielded from light with foil. Avoid exposure to bright, white light. This is because the free chromophore(once**

released from its protein-bound state) is very susceptible to photobleaching. The aim is to exclude/minimize photodamage.

Day 2 - Methanolysis

6 - Resuspend and pool pellets in a final volume of 500 ml MeOH.

7 - Heat suspension in a 1-liter flask in a water bath at 70 - 75 °C with a condensing coil cooled with tap water (see below) for 5-8 hrs (very slow boil).

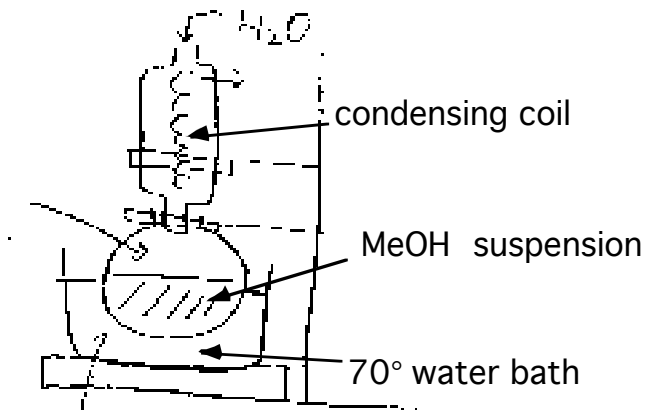

8 - Transfer suspension to two 250 ml bottles and centrifuge 8000 RPM, GS3 rotor, 4° C / 20 min.


9 - Decant and filter supernatant through miracloth directly into a 1-liter side-armed flask and store, foil-wrapped at -20 °C (1st methanolysis supernatant). Store firm pellet from centrifugation, foil-wrapped at -20 °C.

Day 3

10 - Set up 2nd methanolysis of pellets from step 9 as in steps 6-9. (Alternatively, set up at the end of Day 2 and run methanolysis overnight). (2nd methanolysis supernatant).

Day 4

11 – Remove 1st methanolysis supernatant (step 9) from freezer and connect to another side-armed flask which is attached to a vacuum pump and sitting in a dry ice / EtOH bath (cold trap collection flask). Monitor volume of MeOH in collection flask until it reaches 400 ml. This process will take 9-10 hrs depending on vacuum. Store concentrated PCB/MeOH remaining in the original flask (~50 ml) foil-wrapped, at -20 °C.

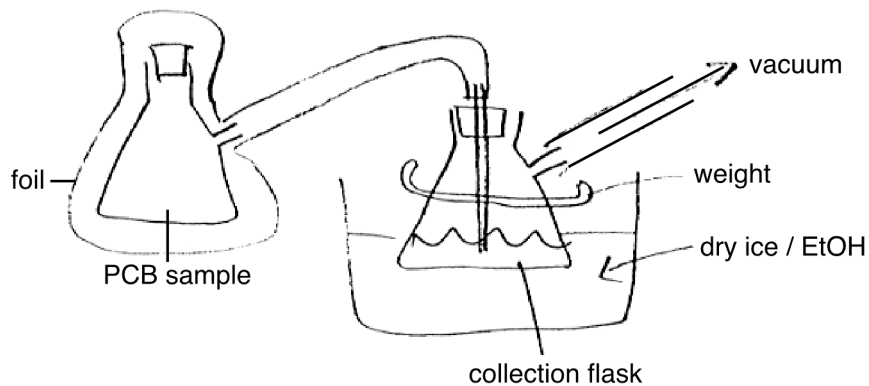

Day 5

Chloroform extraction of phycocyanobilin (PCB)

12 - In a large separatory funnel – mix concentrated PCB/MeOH (~50 ml) with 50 ml chloroform and 100 ml H₂O. Shake and let stand for 5 min.

13 – Collect chloroform (bottom phase) in a beaker. This contains the PCB.

14 - Add another 50 ml chloroform to remaining aqueous phase and repeat steps 13 and 14. Continue with chloroform extraction until solution is almost colorless (usually 2-3X)

15 – Pool these several chloroform/PCB fractions and transfer to a 1 liter round bottom flask. Evaporate chloroform with a stream of nitrogen gas.

16 – Dissolve dried PCB in 4 X 0.75 ml DMSO – Final volume ~3 ml. This is the stock solution. Remove samples to determine PCB concentration

(see below), then aliquot remaining DMSO/PCB and store, foil-wrapped, at -80 °C.

17 – Repeat steps 11-16 on 2nd methanolysis supernatant.

Determination of PCB concentration

Dilute DMSO/PCB stock (1:100, 1:1000) into 1 ml MeOH:HCl, 95%:5% solution for the determination of [PCB] in a spectrophotometer (HCl is 37.5%)


Aliquot DMSO/PCB stock and store, foil-wrapped, at -80° C

Read Absorbance at 680 nm.

$$C(mM) = \frac{A_{680}}{37.9} \times \text{dil. factor}$$

Scan sample 300 – 800 nm for Absorbance Spectrum.

A sample spectrum of a good PCB prep (1:1000) is shown below:

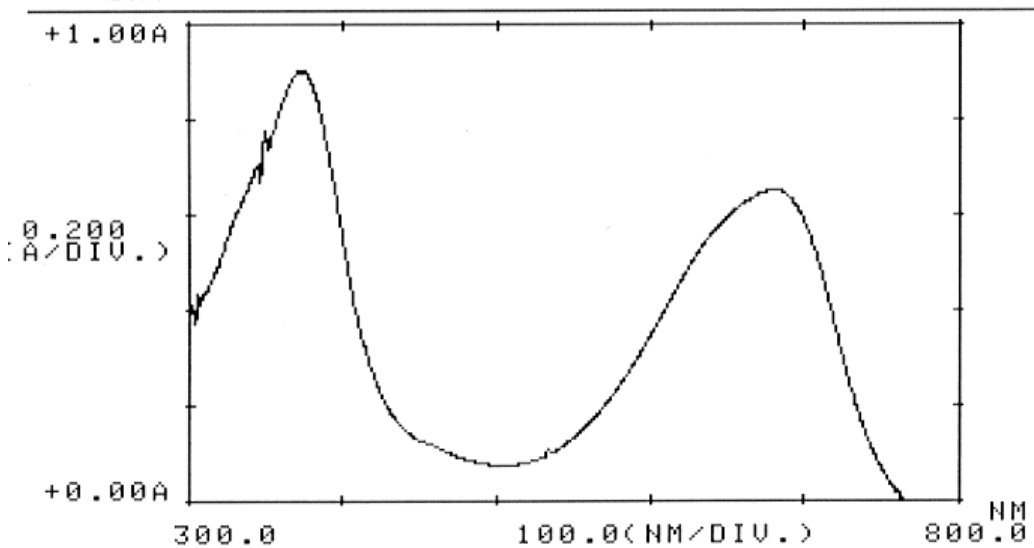

***Safelight Specifications**

Fluorescent tubes: any green fluorescent tube. Example: Sylvania F40G
(www.sylvania.com)

Filters: Roscolene 877 (medium blue green)
Roscolene 874 (medium green)
Thin, flexible colored plastic sheets.
(www.roscow.com)

Green tubes must be wrapped completely with both filters – one layer of each.

Note: This safelight is also used 9in darkroom space for manipulations during experiments with the gene-promoter system in cells.

For questions on this protocol, contact:

Jim Tepperman (Quail Lab)
Plant Gene Expression Center
800 Buchanan St.
Albany, CA 94710


jmtepp@nature.berkeley.edu

510-559-5935


Plasmids for Light-induced Transactivation

PhyA-Full Length and phyB-N-terminal fusions to the GAL4 – binding domain (GBD).


:

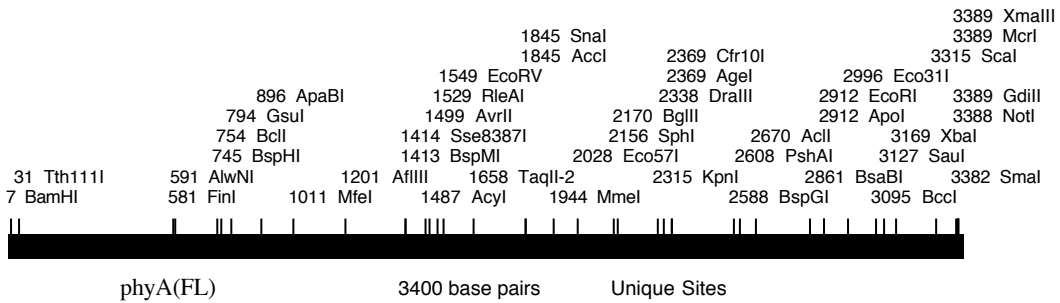

BamHI:NotI fragment in D153 (see below)

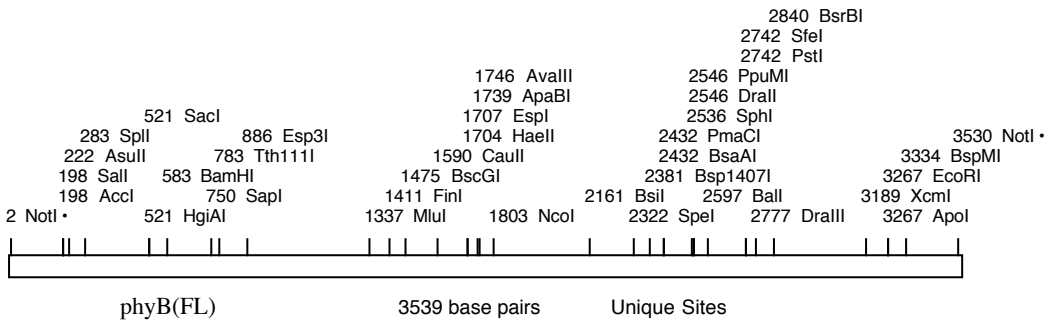

GADPIF3 plasmid:

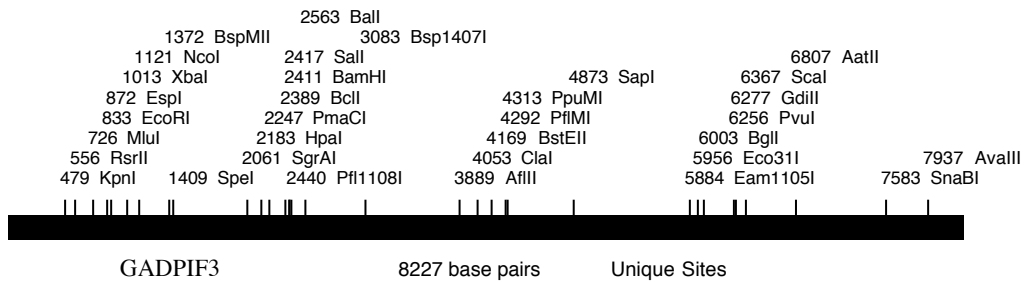

PIF3 ORF: EcoRI (833) – BamHI (2411)

vector; pGAD424 Clontech.

<http://www.clontech.com/techinfo/vectors/vectorsF-I/pGAD424.shtml>